

# DUAL COLOR THC TESTING KIT

Read all SAFETY INFORMATION before use.



**Compassionate  
Analytics**

Plant Knowledge

## SAFETY

- A. These kits use chemicals which may be dangerous to inhale, ingest, or touch. READ ALL MATERIAL SAFETY DATA SHEETS (available for download at [www.compassionateanalytics.com](http://www.compassionateanalytics.com)).
  - ✓ Use in a well ventilated area or a fume hood
  - ✓ Wear gloves at all times
  - ✓ Minimize exposure by keeping lids and caps on tubes as much as possible – only open as necessary, and always hold tubes when open.
- B. RED and YELLOW indicator dyes may stain clothing if spilled – keep paper towel or other absorbent nearby in case of accidental spill.
- C. Disposal (where local regulations allow)
  - ✓ EXTRACTION SOLUTION contains Isopropanol and other nonhazardous ingredients. May be diluted with lots of water and poured down the drain.
  - ✓ RED and YELLOW indicator dyes – after use, dispose of tubes (tightly capped) containing indicator dyes.

## TIPS AND TRICKS

- A. Only 1 Extraction Tube is provided per kit. Rinse and reuse for each test. Contact Compassionate Analytics for additional tubes
- B. For Step 2B – it is important to use a scale that is accurate to 0.01g, as this will affect the accuracy of the test. Contact Compassionate Analytics if you need to obtain an appropriate scale.
- C. During Step 3B – when mixing the ASSAY TUBES, some particulate will remain in the tube. This is normal and will not affect results
- D. The RED and YELLOW indicator dyes may react with some other cannabinoids such as CBD. Most strains have very low levels of other cannabinoids, and thus the results will be unaffected. If you suspect your strain has high CBD, you may use our CBD Test Kit to determine the CBD%.